



SPECIFICATION SHEET

Large Volume UltraMount

INTENDED USE: UltraMount is formulated for covering tissue sections and cell preparations immunohistochemically stained with chromogens producing alcohol soluble OR alcohol insoluble end products.

APPLICATION: UltraMount can be used with chromogens such as Fast-Red or AEC that require an aqueous mounting medium, or those that do not, such as DAB. No heating is required prior to use. This product is optimal for use without cover slips.

AVAILABILITY:

<u>CATALOG #</u>	<u>VOLUME</u>
TA-060-UM	60 ml
TA-125-UM	125 ml

PROTOCOL: Apply several drops of UltraMount to stained specimens. Allow to dry.

STORAGE: Store at room temperature. Do not use past the expiration date.

PRECAUTIONS: Avoid contact with skin and eyes. Wash after use. Wear appropriate protective equipment. Observe all federal, state, and local environmental regulations regarding disposal. Refer to MSDS for additional information.

LIMITATIONS: For Laboratory Use

47790 Westinghouse Dr. Fremont CA 94539 USA
Tel: (800) 828-1628 Fax: (510) 991-2826
Email: labvision@labvision.com Website: www.labvision.com



TA-125-UM
Rev 062703
Page 1 Of 1