



SPECIFICATION SHEET

Fast-Red Substrate System

DESCRIPTION:

Fast-Red in naphthol phosphate is a widely used chromogen for immunohistochemical staining and immunoblotting. In the presence of alkaline phosphatase enzyme, Fast-Red produces a bright red precipitate that is soluble in alcohol.

AVAILABILITY:

<u>CATALOG #</u>	<u>CONTENTS</u>
TA-060-AF	12 Fast Red Tablets 60ml Naphthol Phosphate Substrate
TA-125-AF	25 Fast Red Tablets 125ml Naphthol Phosphate Substrate

IMMUNOHISTOCHEMICAL STAINING PROTOCOL:

1. Add 1 Fast-Red tablet to 5ml of Naphthol Phosphate Substrate and shake until fully dissolved.
2. Apply solution to tissue section.
3. Incubate tissue section for 10-25 minutes.
4. Rinse tissue, counterstain, and coverslip using an aqueous mounting media.

STORAGE:


Store at 2-8°C. Each component is stable for 18 months.

PRECAUTIONS:

47790 Westinghouse Dr. Fremont CA 94539 USA
Tel: (800) 828-1628 Fax: (510) 991-2826
Email: labvision@labvision.com Website: www.labvision.com



TA-060-AF
Rev 21127
Page 1 Of 1

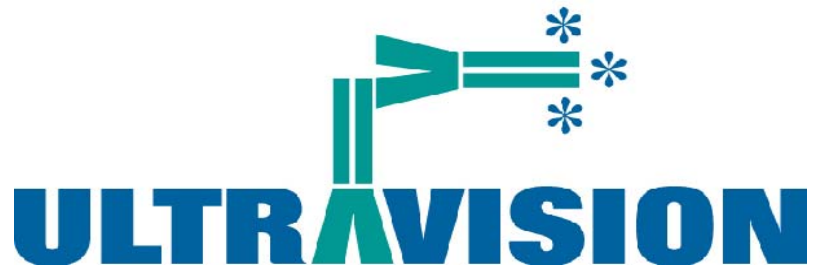

Do not pipet by mouth. Avoid contact with skin and eyes. Handle with care and dispose of according to regulations.

LIMITATIONS:

This product is available for research use only; not for diagnostic or therapeutic work. Lab Vision Corporation will not be held responsible for patent infringement or other violation that may occur with the use of this product.